Instruction Manual
Manuale di istruzioni
Manuel d'instructions
Manual de instrucciones
Bedienungsanleitung
指导手册

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This unit has been designed and manufactured in compliance with the following standards:

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Safety requirements for electrical equipment for measurement, control and for laboratory use

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Electrical equipment for laboratory use

UL 61010-1

General requirement - Canadian electrical code

CAN/CSA-C22.2 No.61010-1

VELP reserves the right to modify the characteristics of its products with the aim to constantly improving their quality.

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It is the responsibility of the user, to properly decontaminate the unit in case of hazardous substances remaining on the surface or interior of the device. Non-reactant decontaminating substances are recommended for use when cleaning, to avoid chemical reaction with the material(s) contained therein. If in doubt about the compatibility of a cleaning or decontamination product, contact the manufacturer or distributor

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The plug disconnects the instrument. Therefore, place the instrument where it can be quickly disconnected.

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Use only power cables provided with three terminals that is wired to ground

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Keep all electrical cords, cables and wires away from liquids.

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Position the instrument on a flat surface, with a distance from the wall of 30 cm (at least).

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If unit is not operating properly do not use. In case of any malfunction, contact your nearest service center

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Do not use with explosive and dangerous materials for which the equipment is not designed. The device must not be used in explosive atmospheres.

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The safety of the instrument is no longer guaranteed in the case of improper use or if the instructions in this manual are not followed.

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JPV is composed by a plastic water tank in which an impeller driven by an electric motor forces the water into four water jets developing suction. The water is continuously recirculated and its consumption is nil.

JPV can be connected to laboratory equipment as rotating evaporators, Buchner, Gooch or membrane filters, gel dryers, etc. The pump produces at the aspiration tubes a residual pressure settable up to 35 mm Hg.

For the movement of equipment use the two handles located in front and rear of the tank. To remove the motor from tank use the two lateral handles located on the plastic cover. In the upper part of the instrument a tube holder connects to the internal Venturi ejectors which produce a maximum air flow of 22 l/minute and a vacuum residual of 35 mm Hg.

The air flow can be regulated by the proper knob. The pump is equipped by a handy door for the introduction of water and a level indicator for maximum and minimum. A stopcock located in the rear part of the tank allows the suggested periodical drainage. On the pipe connecting piece, inside the instrument, there is a safety valve that blocks possible water back-flows in the suction pipe when a pressure differences arise during working cycle (black-out).

2. Assembly and installation

Upon receipt and after having removed the packaging, please check the integrity of the instrument. The box includes:

- JPV recirculating water pump
- Power cable
- Instruction manual

Place the instrument at about 30 cm from the walls.

The equipment is to be connected to the right electric supply by the electric cable connected to the rear of the upper part by a panel socket with fuse. Ensure that the electric mains is earthed and that the mains socket is easy to reach.

Fill the tank with tap water through the front door up to the maximum level shown by the external indicator.

Turn on the instrument by the lighted main switch.

Air flow and vacuum level are regulated by the proper knob.

Air flow: from 15 to 35 l/minute

Vacuum: from room pressure (1 bar) to 0.04 bar

3. Maintenance

In compliance with the product guarantee law, repairs to our units must be carried out in our factory, unless previously agreed otherwise with local distributors. The instrument must be transported in its original packaging and any indications present on the original packaging must be followed (e.g. palletized).

It is the responsibility of the user, to properly decontaminate the unit in case of hazardous substances remaining on the surface or interior of the device. If in doubt about the compatibility of a cleaning or decontamination product, contact the manufacturer or distributor.

- Replace, periodically the tap water in the tank.
 - During operation, if necessary, replace the level of water in the tank up to the maximum level and no more to avoid spillages. A level of water lower than the minimum can compromise pump efficiency.
- During long idle periods it is suggested to empty completely the tank and to wash with tap water.

3.1 Cleaning

Disconnect the unit from the power supply and use a cloth dampened with an non-inflammable non-aggressive detergent.

4. Technical data

Code	F40100541	F40110541
Voltage supply	230V/50-60Hz +/- 10%	115V50/60Hz +/-10%
Fuses	2 x T3,15A AL 250V	2 x T5A AL 250V
Power	370 W	250 W
Dimensions (WxHxD)	250x400x370 mm	
Weight	10 kg (empty tank)	
Flow regulation	from 15 to 35 l/1'	
Vacuum regulation	from 1 to 0.04 bar	
Residual pressure	35 mm Hg with a water temperature in the tank = 15°C	
Liquid volume	Min. 9 L - Max. 11 L	
Enviromental temp. range	from 0 to 40°C	
Max humidity	80%	
Pollution degree EN61010-1	2	
Overvoltage category EN61010-1	II	
Altitude	Up to 2000 m	
Noise	75 dBA	

Equipment intended to be used in wet location

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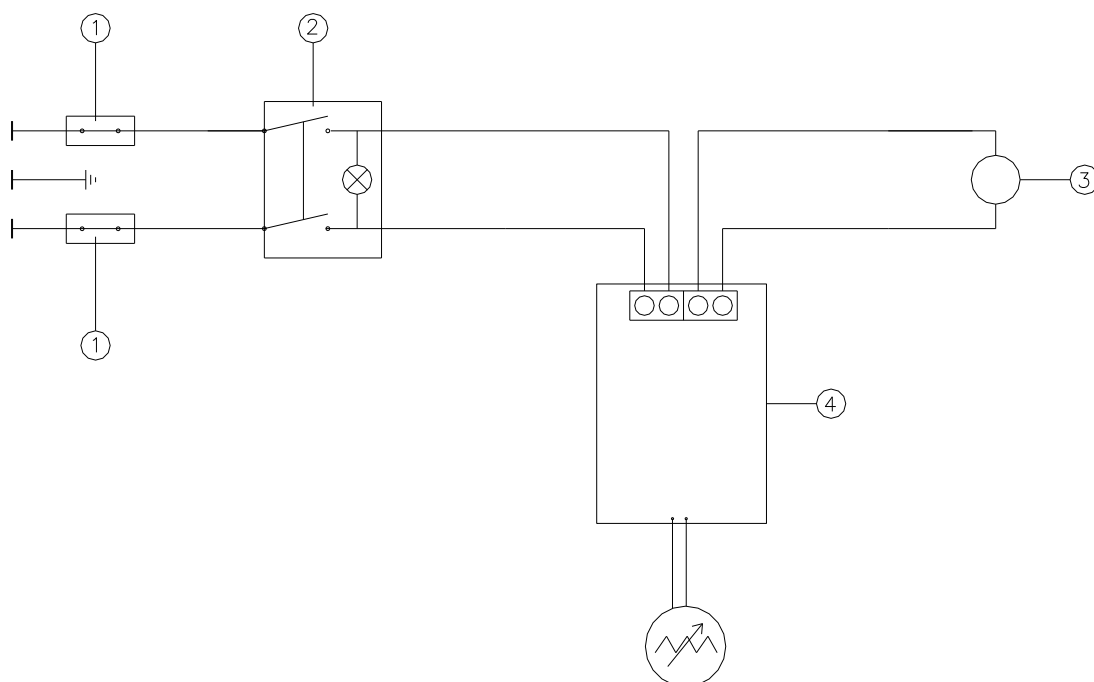
We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following standards:

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EN 61010-1

EN 61326-1

and satisfies the essential requirements of the following directives:

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Machinery directive

2014/30/EU

Electromagnetic compatibility directive

2015/863/EU

(RoHS III)

2012/19/EU

(WEEE)

- plus modifications / più modifiche / plus modifications / más sucesivas modificaciones / in der jeweils gültigen Fassung / 加上修改

7. Declaration of conformity

We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following regulations:

S.I. 2016/1101 Electrical Equipment (Safety) Regulations 2016

S.I. 2016/1091 Electromagnetic Compatibility Regulations 2016

according to the relevant designated standards:

EN 61010-1 Safety requirements for electrical equipment for measurement, control, and laboratory use
Part 1: General requirements

EN 61326-1 Electrical equipment for measurement, control and laboratory use - EMC requirements -
Part 1: General requirements

and satisfies the essential requirements of regulations:

S.I. 2008/1597 Supply of Machinery (Safety) Regulations 2008

S.I. 2012/3032 Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic
Equipment Regulations 2012

S.I. 2013/3113 Waste Electrical and Electronic Equipment Regulations 2013

plus modifications.

Thank you for having chosen VELP!

Established in 1983, VELP is today one of the world's leading manufacturer of analytical instruments and laboratory equipment that has made an impact on the world-wide market with Italian products renowned for innovation, design and premium connectivity. VELP works according to **ISO 9001**, **ISO14001** and **OHSAS 18001** Quality System Certification.

Our instruments are manufactured in Italy according to the IEC 1010-1 and CE regulation.

Our product lines:

Analytical instruments

Elemental Analyzers
Digestion Units
Distillation Units
Solvent Extractors
Fiber Analyzers
Dietary Fiber Analyzers
Oxidation Stability Reactor
Consumables

Laboratory Equipment

Magnetic Stirrers
Heating Magnetic Stirrers
Heating Plates
Overhead stirrers
Vortex mixers
Homogenizers
COD Thermoreactors
BOD and Respirometers
Cooled Incubators
Flocculators
Overhead Shakers
Turbidimeter
Radiation Detector
Open Circulating Baths
Pumps

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VELP Scientifica Srl