

# Lovibond® Water Testing

## Tintometer® Group



### Spectrophotometer XD 7000 (VIS)

Simplify your Work - Save Time and Money!



- Premium optical system with reference beam
- Automatic test recognition with internal barcode reader
- Automatic cuvette type detection
- Support of analytical quality assurance

Part Number: 71307000

#### Latest technology

The XD 7000 and XD 7500 spectrophotometers are equipped with VIS respectively UV / VIS reference beam technology.

#### Automatic test recognition

Each of the more than 150 pre-programmed Lovibond® methods is recognized by barcode as well as the cuvette type.

#### Analytical quality assurance

Standard procedures for analytical quality assurance support the verification of the photometer, the overall system (including the chemical methodology) and the verification of matrix effects.

#### Built-in security levels

The devices support the assignment of passwords and the assignment of up to three different levels of rights.

#### Additional functions

Measurement of transmission and absorption, spectral scan, kinetics analysis, as well as the creation of user-defined methods.

#### Interfaces for data processing

You want to process your data? With Ethernet, USB B, USB A for external memory, keyboard, barcode scanner and printer different versions are available.

#### Industry

Chemical Industry | Food and Beverage Industry | Industries Others | Marine Industry | Municipalities | NGO | Oil and Gas | Pharmaceutical Industry | Power and Energy

#### Application

Boiler Water | Cooling Water | Disinfection Control | Drinking Water Treatment | Food and Beverage | Galvanization | Others | Pool-Water Control | Pool Water Treatment | Raw Water Treatment | Waste Water Treatment

#### Spectrophotometer XD 7000 (VIS)

The XD series instruments are ideal for routine and spectral analysis. As true all-rounders, the XD 7000 and XD 7500 facilitate work routines, especially with automatic method recognition using barcodes, automatic cell recognition and more than 150 pre-programmed methods. In addition to analytical quality control, they also support Good Labor Practice (GLP). The instruments are available together with bar-coded test tubes and a wide range of accessories from a single source - which is also easy on your wallet. High-quality reference beam optics, usability, easy handling and multilingual global applicability round off the versatility of the VIS and UV/VIS spectrophotometers, which are also equipped for flexible on-site use.

## Measuring Range

Test Name	Measuring Range	Chemical Method
Alkalinity-m HR T	5 - 500 mg/L CaCO <sub>3</sub>	Acid / Indicator
Alkalinity-m T	5 - 200 mg/L CaCO <sub>3</sub>	Acid / Indicator
Alkalinity-p T	5 - 500 mg/L CaCO <sub>3</sub>	Acid / Indicator
Aluminium PP	0.01 - 0.25 mg/L Al	Eriochrom Cyanine R
Aluminium T	0.01 - 0.3 mg/L Al	Eriochrom Cyanine R
Ammonia HR TT	1.0 - 50 mg/L N	Salicylate
Ammonia LR TT	0.02 - 2.5 mg/L N	Salicylate
Ammonia PP	0.01 - 0.8 mg/L N	Salicylate
Ammonia T	0.02 - 1 mg/L N	Indophenole Blue
Arsenic	0.02 - 0.6 mg/L As	Silver Diethyldithiocarbamate
Bromine 10 T	0.1 - 3 mg/L Br <sub>2</sub>	DPD
Bromine 50 T	0.05 - 1 mg/L Br <sub>2</sub>	DPD
Bromine PP	0.05 - 4.5 mg/L Br <sub>2</sub>	DPD
Bromine T	0.05 - 13 mg/L Br <sub>2</sub>	DPD
Cadmium M. TT	0.025 - 0.75 mg/L Cd	Cadion
Chloramine (M) PP	0.02 - 4.5 mg/L NH <sub>2</sub> Cl as Cl <sub>2</sub>	Indophenole method
Chloride L (A)	5.00 - 60 mg/L Cl <sup>-</sup>	Iron(III)-thiocyanate
Chloride L (B)	0.5 - 20 mg/L Cl <sup>-</sup>	Mercury Thiocyanate / Iron Nitrate
Chloride T	0.5 - 25 mg/L Cl <sup>-</sup>	Silver Nitrate / Turbidity
Chlorine 10 T	0.1 - 6 mg/L Cl <sub>2</sub>	DPD
Chlorine 50 T	0.02 - 0.5 mg/L Cl <sub>2</sub> <sup>a)</sup>	DPD
Chlorine (free) and Monochloramine	0.02 - 4.50 mg/L Cl <sub>2</sub>	Indophenole method
Chlorine dioxide 50 T	0.05 - 1 mg/l ClO <sub>2</sub>	DPD / Glycine
Chlorine dioxide PP	0.04 - 3.8 mg/l ClO <sub>2</sub>	DPD
Chlorine dioxide T	0.02 - 11 mg/l ClO <sub>2</sub>	DPD / Glycine
Chlorine HR 10 T	0.1 - 10 mg/L Cl <sub>2</sub> <sup>a)</sup>	DPD
Chlorine HR (KI) T (105)	5 - 200 mg/L Cl <sub>2</sub>	KI / Acid
Chlorine L	0.02 - 4.0 mg/L Cl <sub>2</sub> <sup>a)</sup>	DPD
Chlorine MR PP	0.02 - 3.5 mg/L Cl <sub>2</sub> <sup>a)</sup>	DPD
Chlorine PP	0.02 - 2 mg/L Cl <sub>2</sub> <sup>a)</sup>	DPD
Chlorine T	0.01 - 6.0 mg/L Cl <sub>2</sub> <sup>a)</sup>	DPD
Chromium 50 PP	0.005 - 0.5 mg/L Cr <sup>b)</sup>	Diphenylcarbazide
Chromium PP	0.02 - 2 mg/L Cr <sup>b)</sup>	Diphenylcarbazide
COD HR TT	200 - 15000 mg/L COD <sup>b)</sup>	Dichromate / H <sub>2</sub> SO <sub>4</sub>
COD LMR TT	15 - 300 mg/L COD <sup>b)</sup>	Dichromate / H <sub>2</sub> SO <sub>4</sub>
COD LR TT	3 - 150 mg/L COD <sup>b)</sup>	Dichromate / H <sub>2</sub> SO <sub>4</sub>
COD MR TT	20 - 1500 mg/L COD <sup>b)</sup>	Dichromate / H <sub>2</sub> SO <sub>4</sub>
COD VLR TT	2.0 - 60.0 mg/L COD <sup>b)</sup>	Dichromate / H <sub>2</sub> SO <sub>4</sub>
Copper 50 T	0.05 - 1 mg/L Cu <sup>a)</sup>	Biquinoline
Copper L	0.05 - 4 mg/L Cu <sup>a)</sup>	Bicinchoninate
Copper PP	0.05 - 5 mg/L Cu	Bicinchoninate
Copper T	0.05 - 5 mg/L Cu <sup>a)</sup>	Biquinoline
CyA HR T	10 - 200 mg/L CyA	Melamine
Cyanide 50 L	0.005 - 0.2 mg/L CN <sup>-</sup>	Pyridine-barbituric Acid
Cyanide L	0.01 - 0.5 mg/L CN <sup>-</sup>	Pyridine-barbituric Acid
CyA T	10 - 160 mg/L CyA	Melamine
DEHA PP	0.02 - 0.5 mg/L DEHA	PPST
DEHA T (L)	0.02 - 0.5 mg/L DEHA	PPST
Fluoride L	0.05 - 2 mg/L F <sup>-</sup>	SPADNS
Formaldehyde 10 M. L	1.00 - 5.00 mg/L HCHO	H <sub>2</sub> SO <sub>4</sub> / Chromotropic acid
Formaldehyde 50 M. L	0.02 - 1.00 mg/L HCHO	H <sub>2</sub> SO <sub>4</sub> / Chromotropic acid

Test Name	Measuring Range	Chemical Method
Formaldehyde M. TT	0.1 - 5 mg/L HCHO	H <sub>2</sub> SO <sub>4</sub> / Chromotropic acid
H <sub>2</sub> O <sub>2</sub> 50 T	0.01 - 0.5 mg/L H <sub>2</sub> O <sub>2</sub>	DPD / Catalyst
H <sub>2</sub> O <sub>2</sub> HR L	40 - 500 mg/L H <sub>2</sub> O <sub>2</sub>	Titanium Tetrachloride / Acid
H <sub>2</sub> O <sub>2</sub> LR L	1 - 50 mg/L H <sub>2</sub> O <sub>2</sub>	Titanium Tetrachloride / Acid
H <sub>2</sub> O <sub>2</sub> T	0.03 - 3 mg/L H <sub>2</sub> O <sub>2</sub>	DPD / Catalyst
Hardness Ca and Mg L	0.05 - 4 mg/L CaCO <sub>3</sub>	Calmagite
Hardness Ca and Mg MR TT	10 - 360 mg/L CaCO <sub>3</sub>	Calmagite
Hardness Calcium (B) T	20 - 500 mg/L CaCO <sub>3</sub>	Murexide
Hardness Calcium (B) T	50 - 900 mg/L CaCO <sub>3</sub>	Murexide
Hardness total HR T	20 - 500 mg/L CaCO <sub>3</sub> <sup>1)</sup>	Metallphthaleine
Hardness total T	2 - 50 mg/L CaCO <sub>3</sub>	Metallphthaleine
Hazen 24	10 - 500 mg/L Pt	(APHA) Platinum Cobalt Standard Method
Hazen 50	10 - 500 mg/L Pt	(APHA) Platinum Cobalt Standard Method
Hydrazine C	0.01 - 0.7 mg/L N <sub>2</sub> H <sub>4</sub> <sup>c)</sup>	PDMAB
Hydrazine L	5 - 600 µg/L N <sub>2</sub> H <sub>4</sub>	Dimethylaminobenzaldehyde
Hydrazine P	0.05 - 0.5 mg/L N <sub>2</sub> H <sub>4</sub>	Dimethylaminobenzaldehyde
Hypochlorite T	0.2 - 17 % NaOCl	Potassium Iodide
Iron 10 T	0.05 - 1 mg/L Fe	Ferrozine / Thioglycolate
Iron 50 PP	0.01 - 1.5 mg/L Fe <sup>9)</sup>	1,10-Phenanthroline
Iron 50 T	0.01 - 0.5 mg/L Fe	Ferrozine / Thioglycolate
Iron (TPTZ) PP	0.02 - 1.8 mg/L Fe	TPTZ
Iron HR L	0.1 - 10 mg/L Fe	Thioglycolate
Iron in Mo PP (224)	0.01 - 1.8 mg/L Fe	TPTZ
Iron LR L (A)	0.03 - 2 mg/L Fe	Ferrozine / Thioglycolate
Iron LR L (B)	0.03 - 2 mg/L Fe	Ferrozine / Thioglycolate
Iron PP	0.02 - 3 mg/L Fe <sup>9)</sup>	1,10-Phenanthroline
Iron T	0.02 - 1 mg/L Fe	Ferrozine / Thioglycolate
K <sub>S4.3</sub> T	0.1 - 4 mmol/L K <sub>S4.3</sub>	Acid / Indicator
Lead 10	0.1 - 5 mg/L Pb	4-(2-Pyridylazo)-resorcin
Lead (A) TT	0.1 - 5 mg/L Pb	4-(2-Pyridylazo)-resorcin
Lead (B) TT	0.1 - 5 mg/L Pb	4-(2-Pyridylazo)-resorcin
Iodine T	0.05 - 3.6 mg/L I	DPD
Manganese HR PP	0.1 - 18 mg/L Mn	Periodate Oxidation
Manganese L	0.05 - 5 mg/L Mn	Formaloxime
Manganese LR PP	0.01 - 0.7 mg/L Mn	PAN
Manganese T	0.2 - 4 mg/L Mn	Formaloxime
Molybdate HR L	1 - 100 mg/L MoO <sub>4</sub>	Thioglycolate
Molybdate HR PP	0.3 - 40 mg/L Mo	Mercaptoacetic Acid
Molybdate LR PP	0.03 - 3 mg/L Mo	Ternary Complex
Molybdate T	1 - 50 mg/L MoO <sub>4</sub>	Thioglycolate
Nickel 50 L	0.02 - 1 mg/L Ni	Dimethylglyoxime
Nickel L	0.2 - 7 mg/L Ni	Dimethylglyoxime
Nitrate HR	1.2 - 35 mg/L N	2,6-Dimethylphenole
Nitrate LR2 TT	0.2 - 15 mg/L N	2,6-Dimethylphenole
Nitrate LR TT	0.5 - 14 mg/L N	2,6-Dimethylphenole
Nitrate T	0.08 - 1 mg/L N	Zinc Reduction / NED
Nitrate TT	1 - 30 mg/L N	Chromotropic Acid
Nitrite HR PP	2 - 250 mg/L NO <sub>2</sub> <sup>-</sup>	Ferrous Sulfate Method
Nitrite HR TT	0.3 - 3 mg/L N	Sulfanilic / Naphthylamine
Nitrite LR TT	0.03 - 0.6 mg/L N	Sulfanilic / Naphthylamine
Nitrite PP	0.01 - 0.3 mg/L N	Diazotation

Test Name	Measuring Range	Chemical Method
Nitrite T	0.01 - 0.5 mg/L N	N-(1-Naphthyl)-ethylen-diamine
Nitrite VHR L	25 - 2500 mg/L NO <sub>2</sub> <sup>-</sup>	Ferrous Sulfate Method
Oxygen active T	0.1 - 10 mg/L O <sub>2</sub>	DPD
Oxygen dissolved C	10 - 1100 µg/L O <sub>2</sub> <sup>o</sup>	Rhodazine D TM
Ozone 50 T	0.02 - 0.5 mg/L O <sub>3</sub>	DPD / Glycine
Ozone PP	0.015 - 2 mg/L O <sub>3</sub>	DPD / Glycine
Ozone T	0.02 - 2 mg/L O <sub>3</sub>	DPD / Glycine
Phenol T	0.1 - 5 mg/L C <sub>6</sub> H <sub>5</sub> OH	4-Aminoantipyrine
PHMB T	2 - 60 mg/L PHMB	Buffer / Indicator
Phosphate h. TT	0.02 - 1.6 mg/L P <sup>b</sup>	Phosphomolybdenum Blue
Phosphate HR C	1.6 - 13 mg/L P <sup>c</sup>	Vanadomolybdate
Phosphate HR L	5 - 80 mg/L PO <sub>4</sub>	Vanadomolybdate
Phosphate HR T	0.33 - 26.09 mg/L P	Vanadomolybdate
Phosphate HR TT	0.98 - 19.57 mg/L P	Vanadomolybdate
Phosphate LR C	0.016 - 1.6 mg/L P <sup>c</sup>	Stannous Chloride
Phosphate LR L	0.1 - 10 mg/L PO <sub>4</sub>	Phosphomolybic Acid / Ascorbic Acid
Phosphate LR T	0.016 - 1.305 mg/L P	Phosphomolybdenum Blue
Phosphate PP	0.02 - 0.815 mg/L P	Phosphomolybdenum Blue
Phosphate t. TT	0.02 - 1.1 mg/L P <sup>b</sup>	Phosphomolybdenum Blue
Phosphate total HR TT	1.5 - 20 mg/L P <sup>b</sup>	Phosphomolybdenum Blue
Phosphate total LR TT	0.07 - 3 mg/L P <sup>b</sup>	Phosphomolybdenum Blue
Phosphate TT	0.02 - 1.63 mg/L P	Phosphomolybdenum Blue
Phosphonate PP	0.02 - 125 mg/L PO <sub>4</sub>	Persulfate UV Oxidation Method
pH-value HR T	8.0 - 9.6 pH	Thymol Blue
pH value L	6.5 - 8.4 pH	Phenol Red
pH-value LR T	5.2 - 6.8 pH	Bromocresolpurple
pH-value T	6.5 - 8.4 pH	Phenol Red
Polyacrylate L	1 - 30 mg/L Polyacryl	Turbidity
Potassium T	0.7 - 16 mg/L K	Tetraphenylborat Turbidity
SAC 436 nm	0.5 - 50 m <sup>-1</sup>	Direct Reading EN ISO 7887:1994
SAC 525 nm	0.5 - 50 m <sup>-1</sup>	Direct Reading EN ISO 7887:1994
SAC 620 nm	0.5 - 50 m <sup>-1</sup>	Direct Reading EN ISO 7887:1994
Selenium	0.05 - 2 mg/L Se	3,3'-Diaminobenzidine in Toluene
Silcate T	0.05 - 4 mg/L SiO <sub>2</sub>	Silicomolybdenum Blue
Silicate HR PP	1 - 100 mg/L SiO <sub>2</sub>	Silicomolybdate
Silicate L	0.1 - 8 mg/L SiO <sub>2</sub>	Heteropolyblue
Silicate LR PP	0.05 - 1.6 mg/L SiO <sub>2</sub>	Heteropolyblue
Silica VLR PP	0.005 - 0.5 mg/L SiO <sub>2</sub>	Heteropolyblue
Sulphate HR PP	50 - 1000	Bariumsulphate Turbidity
Sulphate PP	5 - 100 mg/L SO <sub>4</sub> <sup>2-</sup>	Bariumsulphate Turbidity
Sulphate T	5 - 100 mg/L SO <sub>4</sub> <sup>2-</sup>	Bariumsulphate Turbidity
Sulphide L	8 - 1400 µg/L S <sup>2-</sup>	Methylene Blue
Sulphide T	0.04 - 0.5 mg/L S <sup>2-</sup>	DPD / Catalyst
Sulphite 10 T	0.1 - 12 mg/L SO <sub>3</sub>	DTNB
Sulphite T	0.1 - 6 mg/L SO <sub>3</sub>	DTNB
Surfactants M. (anion.) TT	0.05 - 2 mg/L SDSA	Methylene Blue

Test Name	Measuring Range	Chemical Method
Surfactants M. (cation.) TT	0.05 - 1.5 mg/L CTAB	Disulphine Blue
Surfactants M. (not ionic) TT	0.1 - 7.5 mg/L Triton X-100	TBPE
Suspended solids 24	10 - 750 mg/L TSS	Turbidity / Attenuated Radiation Method
Suspended solids 50	10 - 750 mg/L TSS	Turbidity / Attenuated Radiation Method
Tannin L	0.5 - 20 mg/L Tannin	
TN HR 2 TT	5 - 140 mg/L N <sup>b</sup> )	2,6-Dimethylphenole
TN HR TT	5 - 150 mg/L N <sup>b</sup> )	Persulphate Digestion
TN LR 2 TT	0.5 - 14 mg/L N <sup>b</sup> )	2,6-Dimethylphenole
TN LR TT	0.5 - 25 mg/L N <sup>b</sup> )	Persulphate Digestion
TOC HR M. TT	50 - 800 mg/L TOC <sup>b</sup> )	H <sub>2</sub> SO <sub>4</sub> / Persulphate / Indicator
TOC LR M. TT	5 - 80 mg/L TOC <sup>b</sup> )	H <sub>2</sub> SO <sub>4</sub> / Persulphate / Indicator
Triazole PP	1 - 16 mg/L Benzotriazole or Tolyltriazole	Catalyzed UV Digestion
Turbidity 50	5 - 500 FAU	Attenuated Radiation Method
Turbidity 24	10 - 1000 FAU	Attenuated Radiation Method
Urea T	0.1 - 2.5 mg/L Urea	Indophenol / Urease
Zinc L	0.1 - 2.5 mg/L Zn	Zincon / EDTA
Zinc T	0.02 - 1 mg/L Zn	Zincon

## Technical Data

<b>Optics</b>	Grid monochromator with reference beam and beam splitter after exit slit
<b>Measurement</b>	Concentration, single and multi-wavelength measurement of absorbance and % transmission, kinetics, spectra
<b>Wavelength Range</b>	320 - 1100 nm
<b>Wavelength Resolution</b>	1 nm
<b>Wavelength Accuracy</b>	± 1 nm on all Holmium peaks
<b>Wavelength Reproducibility</b>	better than 0,5 nm
<b>Spectral Scope</b>	4 nm
<b>Photometric Range</b>	-3.3 - +3.3 Abs
<b>Photometric Resolution</b>	Absorption: 0.001 ; Transmission: 0.1 %
<b>Photometric Accuracy</b>	0.003 Abs below 0.6 Abs ; 0.5 % from 0.6 to 2.0 Abs
<b>Photometric Reproducibility</b>	0.003 Abs below 0.6 Abs ; 0.5 % from 0.6 to 2.0 Abs
<b>Photometric Linearity</b>	< 1 % up to 2.0 Abs between 340 to 900 nm
<b>Scan Speed</b>	700 - 2000 nm/min.
<b>Drift</b>	< 0.005 Abs per hour after 15 minutes heat up time
<b>Stray Light</b>	< 0.1 % Transmission at 340 and 408 nm
<b>Suitable Vials</b>	Rectangular Cuvettes 10 mm Rectangular Cuvettes 20 mm Rectangular Cuvettes 50 mm Round Cuvettes 13 mm Round Cuvettes 16 mm Round Cuvettes 24 mm
<b>Display</b>	7" high contrast colour graphic-display
<b>Interfaces</b>	Ethernet USB B USB A for External Memory Keypad Barcode-Scanner PCL Compatible Printer
<b>Operation</b>	Membrane Keyboard
<b>Automatic Cuvette Recognition</b>	Round cuvettes: 13, 16 and 24 mm ; Rectangular cuvettes: 10, 20 and 50mm

<b>Auto – OFF</b>	Yes
<b>Test Recognition</b>	via internal barcode reader
<b>Auto-Check</b>	Self-test at each switch-on: test of memory, processor, internal interface, filter lamp and additional calibration of each wavelength
<b>LIMS Compatibility</b>	ASCII, .csv-files
<b>Internal Storage</b>	approx. 5000 data sets (method, user ID, date, result), autostorage function / manual storage function
<b>Security</b>	Password protection possible: 3 different user levels (guest, user, admin)
<b>Power Drain</b>	100 - 240 V, 50/60 Hz
<b>Power Supply</b>	Buffer batteries (4 x AA), power supply unit with cable
<b>Portability</b>	Benchtop
<b>Environmental Conditions</b>	+10 °C to 35 °C (41 °F to 95 °F), ≤ 75 % average humidity within one year 95 %, max. 30 days/year, 85 % all other days
<b>Stock Conditions</b>	-25 °C to +65 °C (-13 °F to 268 °F)
<b>Protection Class</b>	IP 30
<b>Compliance</b>	CE
<b>IP Protection Class</b>	EN 60529
<b>Interference Emission</b>	Class B
<b>Interference Immunity</b>	IEC 61000-4-3
<b>Tolerance Extension</b>	0.008 E
<b>Meter Safety</b>	EC Directive 2014/35/EC EN 61010-1:2010
<b>Languages User Interface</b>	German, English, French, Spanish, Italian, Portuguese, Polish, Indonesian, Russian, Chinese, Japanese, Dutch, Swedish, Norwegian, Czech, Romanian, Macedonian, Slovenian, Hungarian, Turkish, Korean, Vietnamese, Thai, Serbian, Malaysian, Danish, Bulgarian
<b>Languages Quick Start Guide</b>	German, English, French, Spanish, Italian, Portuguese, Polish, Indonesian, Russian, Chinese, Japanese, Dutch, Swedish, Norwegian, Czech, Romanian, Macedonian, Slovenian, Hungarian, Turkish, Korean, Vietnamese, Thai, Serbian, Malaysian, Danish, Bulgarian
<b>Languages Full User Manual</b>	German, English, Spanish, French, Italian, Portuguese, Chinese, Japanese
<b>Dimensions</b>	422 x 195 x 323 mm

## Delivery Scope

- 4 batteries (AA)
- 1 power supply cable
- 4 round cuvettes with lid and 1 zero cuvette XD 7x00 (ø 24 mm)
- 1 zero cuvette (ø 16 mm)
- Quickstart guide in 24 languages
- Full user manual in 8 languages
- Warranty information

## Accessories

Title	Part Number
Adapter (13 mm) MultiDirect for Vacu-vial	192075
Batteries (AA), set of 4	1950025
Round cuvette 24 mm, set of 12	197620
Round cuvette 24 mm, set of 5	197629
Cleaning cloth	197635
Round cuvette 16 mm, set of 10	197665
Adapter for round cuvettes 13 mm	19802192
Mixing cylinder, 25 ml	19802650
Inspection and calibration package for XD7000	19802707
Zero cuvette ø 16 mm for XD 7000/7500	215661
Zero cuvette ø 24 mm for XD 7000/7500	215662
Thermoreactor RD 125	2418940
USB cable 3 m	2444482
Pipette, 1000 µl	365045
Measuring spoon, 1 g	384930
UV Pen Lamp, 254 nm	400740
UV protection glasses, orange	400755
Cuvette stand for 6 round cuvettes Ø 24 mm	418951
Cuvette stand for 10 round cuvettes Ø 16 mm	418957
Pipette tips, 1-5 ml (white) 100 pc.	419066
Pipette tips, 0,1-1 ml (white), 1000 pc.	419073
Automatic pipette, 1-5 ml	419076
Automatic pipette, 0,1-1 ml	419077
Screw caps TOC	420757
Measuring spoon no. 8, black	424513
Universal Container - Cap	424648
Plastic funnel with handle (white)	471007
ValidCheck Chlorine 1,5 mg/l	48105510
Stirring rod and spoon	56A006601
W100/OG/10MM Rectangular cell, optical glass	601040
W100/OG/20MM Rectangular cell, optical glass for determination of arsenic	601050
W100/OG/50MM Rectangular cell, optical glass	601070
W110/UV/10MM Rectangular cell, Quartz UV	661130
W110/UV/20MM Rectangular cell, Quartz UV	661140
W110/UV/50MM Rectangular cell, Quartz UV	661160
Energy station XD series/SpectroDirect	711051
Secondary standard set VIS with DAKkS calibration certificate	711160
Light source replacement XD 7000	71310000
Carrying case for XD 7000/7500	71310010
12 V-plug connector for XD 7000/7500	71310020
Barcode Hand-held Scanner	71310030
Semimicro cell, 50 mm with lid	71310045
Factory calibration certificate ISO 9001 for XD7000	999754

---

**Tintometer GmbH**

Lovibond® Water Testing  
Schleefstraße 8-12  
44287 Dortmund  
Tel.: +49 (0)231/94510-0  
sales@lovibond.com  
www.lovibond.com  
Germany

**The Tintometer Limited**

Lovibond House  
Sun Rise Way  
Amesbury, SP4 7GR  
Tel.: +44 (0)1980 664800  
Fax: +44 (0)1980 625412  
sales@lovibond.uk  
www.lovibond.com  
UK

**Tintometer China**

Room 1001, China Life Tower  
16 Chaoyangmenwai Avenue,  
Beijing, 100020  
Tel.: +86 10 85251111 App. 330  
Fax: +86 10 85251001  
chinaoffice@tintometer.com  
www.lovibond.com  
China

**Tintometer South East Asia**

Unit B-3-12, BBT One Boulevard,  
Lebuh Nilam 2, Bandar Bukit Tinggi,  
Klang, 41200, Selangor D.E  
Tel.: +60 (0)3 3325 2285/6  
Fax: +60 (0)3 3325 2287  
lovibond.asia@tintometer.com  
www.lovibond.com  
Malaysia

**Tintometer Brazil**

Caixa Postal: 271  
CEP: 13201-970  
Jundiaí – SP  
Tel.: +55 (11) 3230-6410  
sales@tintometer.com.br  
www.lovibond.com.br  
Brazil

**Tintometer Inc.**

6456 Parkland Drive  
Sarasota, FL 34243  
Tel: 941.756.6410  
Fax: 941.727.9654  
sales@lovibond.us  
www.lovibond.com  
USA

**Tintometer India Pvt. Ltd.**

Door No: 7-2-C-14, 2<sup>nd</sup>, 3<sup>rd</sup> & 4<sup>th</sup> Floor  
Sanathnagar Industrial Estate,  
Hyderabad, 500018  
Telangana  
Tel: +91 (0) 40 23883300  
Toll Free: 1 800 599 3891/ 3892  
indiaoffice@lovibond.in  
www.lovibondwater.in  
India

**Tintometer Spain**

Postbox: 24047  
08080 Barcelona  
Tel.: +34 661 606 770  
sales@tintometer.es  
www.lovibond.com  
Spain

Technical changes without notice  
Printed in Germany

Lovibond® and Tintometer® are Trademarks of the Tintometer Group of Companies