

RapidChek® SELECT

Salmonella Group B

RapidChek® SELECT™ Salmonella Group B test system detects Salmonella **Group B serovars such as** S. Typhimurium and S. Heidelberg in poultry house environmental samples in a fast and easy to use way. The method is a useful serogroup screening tool for poultry processors.



How the Test Works

The RapidChek® technology couples a sensitive immune-detection strip for Salmonella Group B serovars with the SELECT™ Salmonella enrichment media for fast and accurate results within 32 - 44 hours for poultry house environmental samples. The method uses a novel application of bacteriophages in the media to act as selective agents, enhancing both specificity and sensitivity of the overall method. The immune-detection strip employs proprietary anti-Salmonella B serogroup antibodies. The immunodetection strip is added directly into the provided tube with the secondary enrichment sample. Within 10 minutes results are read. One line indicates a negative result and two lines indicate a presumptive positive result. The test kits are stored at room temperature.

Applications

RapidChek® SELECT™ Salmonella Group B method is designed to detect Group B Salmonella species in poultry house environmental samples. The test kit permits the presumptive detection and identification of the target serogroup in 32 - 44 hours. The test can be used in conjuction with the RapidChek® SELECT™ Salmonella species test and/or Salmonella Enteritidis test method or as a standalone method for serogrouping results.

Confirmation

Presumptive positive results must be confirmed by a cultural reference method (USDA MLG, FDA BAM or ISO). At least two different types of selective agars should be plated for best results. RapidChek® SELECT™ Salmonella secondary media samples used in the test procedure can be used for confirmation.

CUSTOMER BENEFITS

Fast & Simple Procedure

- Convenient serogroup screening information
- Days faster than traditional serology
- Minimal training
- · No additional equipment

Easy Resource Management


- · Unlimited scalability
- Storage at room temperature
- · Long shelf life




RapidChek® SELECT™ Salmonella Group B

Pathogen Screening Quick Reference

Read the Package Insert instructions completely before performing any test.

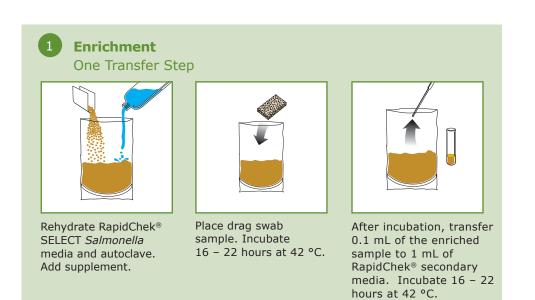

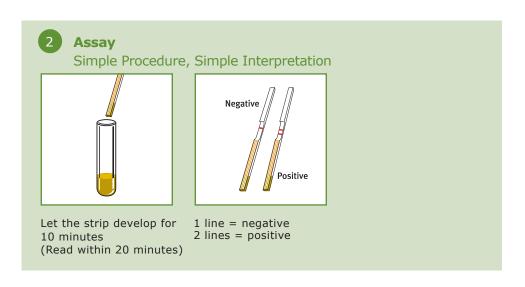

Ordering Information-RapidChek® SELECT™ Salmonella Group B

Item	Description	Part Number
Test Kit	50 Test strips with tubes and pipettes, Media not included	10001419
Media System	500g Primary Enrichment Media, Supplement, 10 g Secondary Media	10001384

Also available:

RapidChek® *E. coli* O157, RapidChek® SELECT™ *Salmonella*, RapidChek® SELECT™ *Salmonella* Enteritidis, RapidChek® *Listeria*, RapidChek® *Listeria* NextDay™, RapidChek® CONFIRM™ non-O157 STEC IMS RapidChek® CONFIRM™ Salmonella Enteritidis IMS

Romer Labs Inc. - America · Newark, DE, USA · T +1 302 781 6400 · F +1 302 781 6378 · E office@romerlabs.com
Romer Labs Diagnostic GmbH - Europe · Tulln, Austria · T +43 2272 615 33 · F +43 2272 615 33 13177 · E office-europe@romerlabs.com
Romer Labs Singapore Pte. Ltd. - Asia · Singapore · T +65 6631 8018 · F +65 6275 5584 · E salesasia@romerlabs.com