

RapidChek® *Campylobacter* Quick Guide

RapidChek® *Campylobacter* Test Kit, 10005433

Test Kit Includes:

- 50 test strips in a desiccated canister
- 50 transfer pipettes
- 50 sample tubes
- Package Insert

Does not include media.

RapidChek® *Campylobacter* Media

- 500 g – 10006043
- 5 kg – 10006042

RapidChek® *Campylobacter* Strips, 10006045

Test Kit Includes:

- 10 x 50 test strips in desiccated canisters

Does not include tubes, pipettes, or media.

Also Available:

- RapidChek® SELECT™ *Salmonella*
- RapidChek® SELECT™ *Salmonella* Enteritidis
- RapidChek® CONFIRM™ *Salmonella* Enteritidis IMS
- RapidChek® *Listeria*
- RapidChek® *Listeria* NextDay™
- RapidChek® *E. coli* O157 (including H7)
- RapidChek® CONFIRM™ non-O157 STEC IMS

RapidChek® *Campylobacter* Comb System, 10006041

Test Kit Includes:


- 8 desiccated foil bags containing 6 combs of 8-strips each
- 48 strips of cluster tubes, 8 tubes x 1.1 mL
- Cluster tube rack, 96-position
- Package Insert

Does not include pipettes or media.



RapidChek®. Simply Accurate.

RapidChek® *Campylobacter* Quick Guide

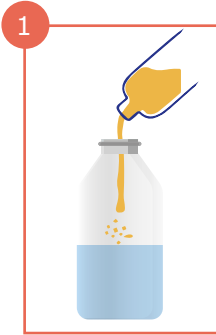

Media Preparation

Add media to sterile water. DO NOT AUTOCLAVE.

Sterile Water (Liters)	RapidChek® <i>Campylobacter</i> Media (Grams)
1	55.2

Rehydrated Media Shelf-Life


6 hours at room temperature.



Sample Enrichment

Add prepared RapidChek® *Campylobacter* Media to the carcass rinse. Incubate at 42 °C in a vented culture flask.

Sample Type	Volume of Media to add (mL)	Incubation Time (Hours)
30 mL carcass rinse	30 mL	47 - 49

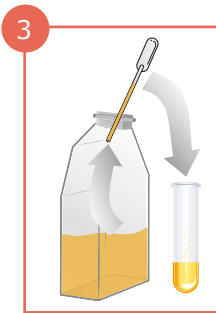


Transfer

Transfer 0.4 mL of the enriched sample into a supplied tube.



Test

Place strips into tubes with arrows facing down.

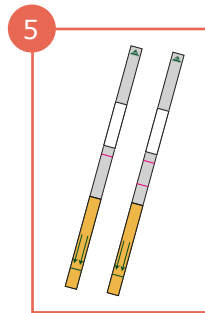

Read Result

After 20 minutes, read result.
1 line = Negative
2 lines = Positive