


# Lovibond® Colour Measurement Tintometer® Group

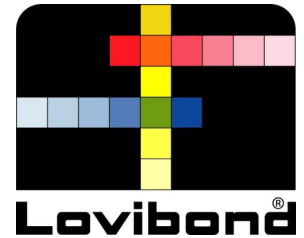

## Model F

Visual Colorimetry : Lovibond® Precision



- Glass Colour Standards Housed in Durable Racks
- Prismatic Optical System Provides Improved Colour Matching
- Balanced Field of View
- Replaceable Sample Chamber Liner

Part Number: 180000

### Visual Colorimeter for Lovibond® RYBN Scale

The Model F is a visual colorimeter designed to optimise the use of Lovibond® glass filters in order to express colour in terms of Lovibond® Red, Yellow, Blue and Neutral.

### Two Adjacent Fields of View

It is arranged with two adjacent fields of view, seen through the viewing tube, so that the product in the sample field and a white reflective surface in the comparison field are observed side by side, suitably illuminated.

### Easy Colour Comparison

The Lovibond® glasses are introduced into the comparison field by a simple system of sliding racks, allowing you to compare the colour of light which is either transmitted through, or reflected from, the sample. The racks are varied until a visual colour match is found for the light from the sample and its colour can then be expressed in Lovibond® units.

### Excellent Colour Matching

A series of neutral glasses in racks is also supplied; these can be introduced into the sample field to 'dull' the colour of products which are too bright to obtain a good colour match using Lovibond® Red, Yellow or Blue glasses. The racks allow a visual colour match to be found from the light from the sample and its colour can then be expressed in Lovibond® units.

### Assured Accuracy of Lovibond® RYBN Colour

Lovibond® Tintometer® is the only company that can guarantee accurate, validated Lovibond® RYBN Colour readings.

### Easy Maintenance for Optimum Performance

The Lovibond® Tintometer is a precision optical instrument which should be kept clean to preserve its accuracy, even in the harsh operating conditions often encountered. This is why the Model F has been designed to allow easy cleaning and maintenance with the minimum of cost and effort. The instrument includes a white plastic liner in the sample chamber which can be removed for cleaning if a spillage takes place. It can be replaced at intervals to maintain the whiteness of the interior, an essential feature for accurate colour matching.

Separate filters in the racks also simplify cleaning and allow replacement of single glasses (as opposed to complete racks) if required.

### Application Versatility

The Model F is an easy-to-use instrument with versatile application in colour measurement of products which transmit light as well as opaque solids, powders and pastes (an optional 'solid' sample accessory pack may be required for colour grading of light-reflecting products).

### Compliance to International Standards

It is available in slightly varying formats to meet the requirements of different applications and the national and international standardising bodies which specify the instrument in their official methods for colour measurement.

## Industry

Chemical Industry | Food and Beverage Industry | Industries Others | Oil and Gas

## Application

Beer, Malts and Caramels | Edible Oils and Fats | Food and Beverage | Petroleum Oils and Waxes

## Model F

The Model F is a visual colorimeter designed to optimise the use of Lovibond® glass filters in order to express colour in terms of Lovibond® Red, Yellow, Blue and Neutral. It is available in slightly varying formats to meet the requirements of different applications and the national and international standardising bodies which specify the instrument in their official methods for colour measurement.

## Technical Data

<b>Applicable Standard</b>	AOCS Cc 13e-92 AOCS Cc 13j-97 BS684 IS 548 (part 1)-2010 ISO 15305 ISO 27608 MS 252 : Part 16 MS 817 : Part 12
<b>Color Type</b>	Transparent Translucent Opaque
<b>Path Length</b>	Up to 153 mm (6")
<b>Illuminants</b>	Approximates to daylight
<b>Portability</b>	Benchtop
<b>Compliance</b>	CE
<b>Dimensions</b>	330 x 230 x 410 mm

## Delivery Scope

- Model F Housing
- Complete set of 11 racks containing permanently coloured glass filters for the Lovibond® colour scale
  - (Red 0.1 - 0.9, 1.0 - 9.0, 10.0 - 70; Yellow 0.1 - 0.9, 1.0 - 9.0, 10.0 - 70; Blue 0.1 - 0.9, 1.0 - 9.0, 10.0 - 40; Neutral 0.1 - 0.9, 1.0 - 3.0)
- Sample chamber liner with a white PTFE powder reference
- Spare white reference
- A pair of spare bulbs
- 1 x 1" & 1 x 5¼" rectangular fused glass cells

## Accessories

Title	Part Number
Lamp 12v10w	124100
Vacuum formed tray	180110
Halon Self Adhesive	183079

**Tintometer GmbH**  
Lovibond® Water Testing  
Schleefstraße 8-12  
44287 Dortmund  
Tel.: +49 (0)231/94510-0  
sales@lovibond.com  
www.lovibond.com  
Germany

**The Tintometer Limited**  
Lovibond House  
Sun Rise Way  
Amesbury, SP4 7GR  
Tel.: +44 (0)1980 664800  
Fax: +44 (0)1980 625412  
sales@lovibond.uk  
www.lovibond.com  
UK

**Tintometer China**  
Room 1001, China Life Tower  
16 Chaoyangmenwai Avenue,  
Beijing, 100020  
Tel.: +86 10 85251111 App. 330  
Fax: +86 10 85251001  
chinaoffice@tintometer.com  
www.lovibond.com  
China

**Tintometer South East Asia**  
Unit B-3-12, BBT One Boulevard,  
Lebuh Nilam 2, Bandar Bukit Tinggi,  
Klang, 41200, Selangor D.E  
Tel.: +60 (0)3 3325 2285/6  
Fax: +60 (0)3 3325 2287  
lovibond.asia@tintometer.com  
www.lovibond.com  
Malaysia

**Tintometer Brazil**  
Caixa Postal: 271  
CEP: 13201-970  
Jundiaí – SP  
Tel.: +55 (11) 3230-6410  
sales@tintometer.com.br  
www.lovibond.com.br  
Brazil

**Tintometer Inc.**  
6456 Parkland Drive  
Sarasota, FL 34243  
Tel: 941.756.6410  
Fax: 941.727.9654  
sales@lovibond.us  
www.lovibond.com  
USA

**Tintometer India Pvt. Ltd.**  
Door No: 7-2-C-14, 2<sup>nd</sup>, 3<sup>rd</sup> & 4<sup>th</sup> Floor  
Sanathnagar Industrial Estate,  
Hyderabad, 500018  
Telangana  
Tel: +91 (0) 40 23883300  
Toll Free: 1 800 599 3891/ 3892  
indiaoffice@lovibond.in  
www.lovibondwater.in  
India

**Tintometer Spain**  
Postbox: 24047  
08080 Barcelona  
Tel.: +34 661 606 770  
sales@tintometer.es  
www.lovibond.com  
Spain

Technical changes without notice  
Printed in Germany

Lovibond® and Tintometer® are Trademarks of the Tintometer Group of Companies